


# ATR – Theory and Applications

Attenuated Total Reflectance (ATR) is today the most widely used FTIR sampling tool. ATR generally allows qualitative or quantitative analysis of samples with little or no sample preparation, which greatly speeds sample analysis. The main benefit of ATR sampling comes from the very thin sampling pathlength and depth of penetration of the IR beam into the sample. This is in contrast to traditional FTIR sampling by transmission where the sample must be diluted with IR transparent salt, pressed into a pellet or pressed to a thin film, prior to analysis to prevent totally absorbing bands in the infrared spectrum.


A comparison of transmission versus ATR sampling results for a thick polymer sample is shown below where the sample is too thick for high-quality transmission analysis (shown in the lower blue spectrum). In transmission spectroscopy, the IR beam passes through the sample and the effective pathlength is determined by the thickness of the sample and its orientation to the directional plane of the IR beam. Clearly in the example below the sample is too thick for transmission analysis because most of the IR bands are totally absorbing. However, simply placing the thick sample on the ATR crystal (Diamond MIRacle) and applying pressure generates a nearly perfect result (upper red spectrum) – identified by a library search as a polybutylene terephthalate. The total analysis time for the thick polymer by ATR was less than 1 minute.

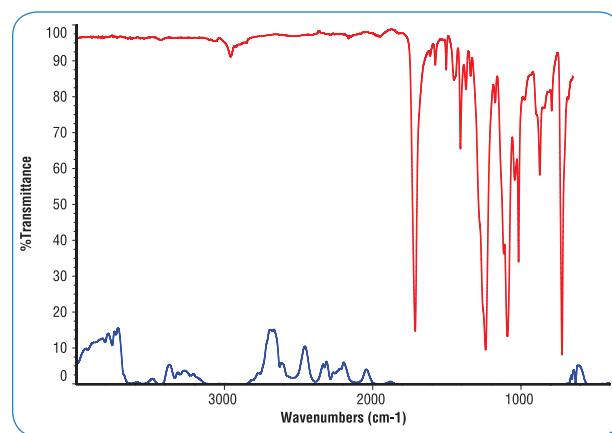

ATR and transmission spectra of a thick polymer sample.

## How ATR Works

With ATR sampling we direct the IR beam into a crystal of relatively higher refractive index. The IR beam reflects from the internal surface of the crystal and creates an evanescent wave, which projects orthogonally into the sample in intimate contact with the ATR crystal. Some of the energy of the evanescent wave is absorbed by the sample and the reflected radiation is returned to the detector. This ATR phenomenon is shown graphically in the following representation of a single reflection ATR.

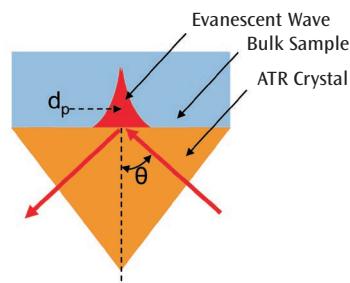

Graphical representation of a single reflection ATR.

While the analysis of samples by ATR is easy, it is interesting and useful to be aware of each of the following experimental factors and how they affect the final spectrum:


- Refractive indices of the ATR crystal and the sample
- Angle of incidence of the IR beam
- Critical angle
- Depth of penetration
- Wavelength of the IR beam
- Effective pathlength
- Number of reflections
- Quality of the sample contact with ATR crystal
- ATR crystal characteristics

The **refractive indices of the crystal and sample** are important considerations in the ATR sampling technique by virtue of the following equation:

$$\theta_c = \sin^{-1} \left( \frac{n_2}{n_1} \right)$$

where  $\theta_c$  is the **critical angle**,  $n_2$  is the refractive index of the sample and  $n_1$  is the refractive index of the crystal.

When the angle of incidence exceeds the critical angle, we will observe a purely ATR spectral result. If the critical angle is not met, we will observe a combined ATR and external reflectance result. This occurs if the angle of incidence of the IR beam is too low, if the refractive index of the crystal is too low, if the refractive index of the sample is too high or a combination of these three factors. In most cases this problem is not observed; however, an example of this is shown in the following spectral data. The sample is a high refractive index liquid ( $n_s=1.8$ ) run on a 45 degree accessory using diamond and Ge crystal plates. The spectrum run on the Ge crystal plate exhibits a normal baseline and symmetric absorbance bands – critical angle is met. The spectrum run on the diamond crystal plate has a baseline shifted and asymmetric absorbance bands due to non-adherence to the critical angle requirements for this set of analysis parameters.

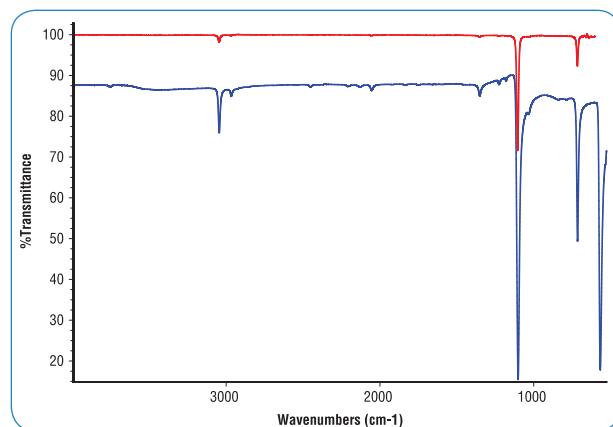

Spectra of high refractive index liquid using Ge (red) and diamond (blue) ATR crystals.

Another way to correct the spectral artifacts observed (above) in the high refractive index sample spectrum would be to increase the angle of incidence in the ATR accessory to a value above the critical angle. Adjustment or selection of the angle of incidence is available in several of the PIKE Technologies ATR accessories.

**Table 1: Pathlengths and penetration depth (in microns), and ATR crystals at various angles of incidence (in degrees) and numbers of reflections.**

ATR Sampling for $n_2 = 1.5$ $\lambda = 1000 \text{ cm}^{-1}$		ZnSe, Diamond $n_1 = 2.4$ $\theta_c = 38.7^\circ$			AMTIR $n_1 = 2.5$ $\theta_c = 36.9^\circ$			Ge $n_1 = 4.0$ $\theta_c = 22.0^\circ$		
$\theta$	$N$	$d_p$	$d_e$	EPL	$d_p$	$d_e$	EPL	$d_p$	$d_e$	EPL
45	1	2.0	4.36	4.36	1.7	3.38	3.38	0.66	0.61	0.61
45	3	2.0	4.36	13.08	1.7	3.38	10.15	0.66	0.61	1.84
45	10	2.0	4.36	43.60	1.7	3.38	33.84	0.66	0.61	6.14
30	1	N/A	N/A	N/A	N/A	N/A	N/A	1.2	1.59	1.59
30	3	N/A	N/A	N/A	N/A	N/A	N/A	1.2	1.59	4.76
30	10	N/A	N/A	N/A	N/A	N/A	N/A	1.2	1.59	15.85
60	1	1.11	1.53	1.53	1.02	1.30	1.30	0.51	0.32	0.32
60	3	1.11	1.53	4.59	1.02	1.30	3.91	0.51	0.32	0.97
60	10	1.11	1.53	15.32	1.02	1.30	13.03	0.51	0.32	3.23

Note: N/A indicates critical angle is violated.

Further useful consideration for ATR analysis is the **depth of penetration** ( $d_p$ ) of the IR beam into the sample. Technically, this is defined as the distance required for the electric field amplitude to fall to  $e^{-1}$  of its value at the surface and is further defined by

$$d_p = \frac{\lambda}{2\pi(n_1^2 \sin^2 \theta - n_2^2)^{1/2}}$$

where  $\lambda$  is the wavelength of light and  $\theta$  is the angle of incidence of the IR beam relative to a perpendicular from the surface of the crystal. Typical depth of penetration in ATR ranges from about 0.5 microns up to about 5 microns depending upon these experimental conditions. Shown in the graphical representation of the ATR phenomenon, the strength of the evanescent wave decays rapidly as we progress from the surface of the ATR crystal. If we wish to compare the sample absorbance of the ATR measurement with that of a transmission measurement, we need to calculate the volume of the evanescent wave, known as the effective penetration of the IR beam. The effective penetration ( $d_e$ ) is unique for parallel polarization ( $d_{e\parallel}$ ) and perpendicular polarization ( $d_{e\perp}$ ) and these are defined by:

$$d_{e\perp} = \frac{n_1^2 n_2 \cos \theta}{(n_1^2 - n_2^2)} \times \frac{\lambda_1}{\pi \sqrt{n_1^2 \sin^2 \theta - n_2^2}}$$

$$d_{e\parallel} = \frac{n_1^2 n_2 \cos \theta}{(n_1^2 - n_2^2)} \times \frac{2n_1^2 \sin^2 \theta - n_2^2}{(n_1^2 - n_2^2) \sin^2 \theta - n_2^2} \times \frac{\lambda_1}{\pi \sqrt{n_1^2 \sin^2 \theta - n_2^2}}$$

$$\text{Where } \lambda_1 = \lambda / n_1$$


The effective penetration for an unpolarized IR beam is the average of the parallel and perpendicular penetration.

$$d_e = \frac{(d_{e\perp} + d_{e\parallel})}{2}$$

Generally, a single reflection ATR is ideal for qualitative analysis, "what is my sample?" When we need to look at minor components of a sample for qualitative or quantitative analysis, then we need to increase the **effective pathlength** (EPL) by increasing the number of reflections ( $N$ ) within the ATR crystal. The effective pathlength in ATR is derived by the following equation, where  $N$  = number of reflections on the sample.

$$EPL = N \times d_e$$


An example of the benefit of increased number of reflections is shown in the following spectral data for the analysis of carbohydrate content in a soft drink sample. The red spectrum is run using a 10-reflection HATR accessory. The blue spectrum is run using a single-reflection ATR using an identical scaling factor. Clearly the minor carbohydrate bands are more readily apparent in the multi-reflection ATR accessory.

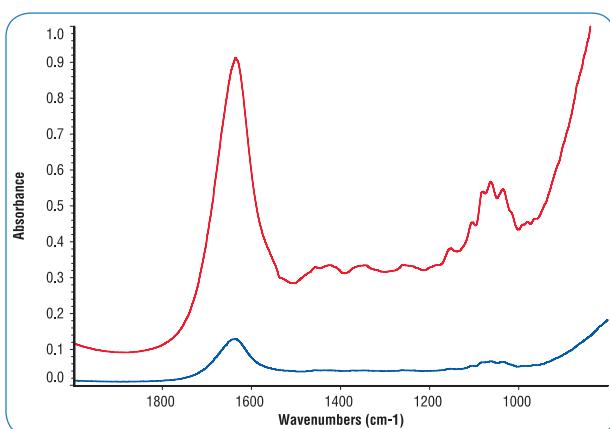

Soft drink sample using 10-reflection and 1-reflection ATR.

For your convenience we have calculated theoretical values of depth of penetration, effective penetration, and effective pathlength for typical combinations of crystal materials, angles of incidence, and number of reflections (Table 1).

With the thin penetration of the evanescent wave into the sample, it is obvious that intimate contact of the sample be made onto the surface of the ATR crystal. For liquid or pliable samples, **quality of sample contact with the ATR crystal** is generally not a problem. For rigid, irregular shaped or porous samples, high pressure sufficient to deform the sample will increase the extent of sample contact and thereby increase sample absorbance. This is shown in the following spectral data collected for a porous foam polymer using a MIRacle ATR with ZnSe crystal.

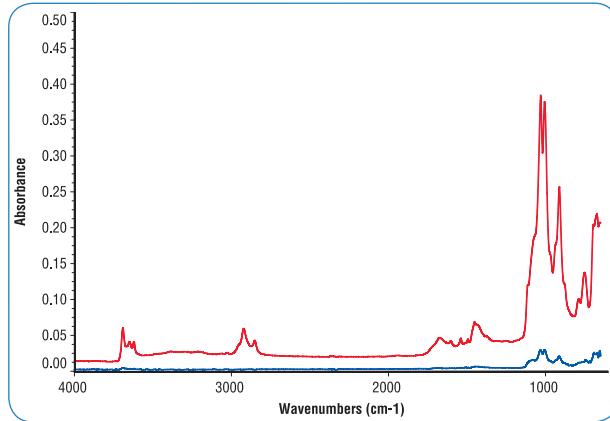

Porous foam sample with high pressure (red) and low pressure (blue).

The blue spectrum was collected with low pressure applied to the foam sample, whereas the red spectrum is produced with high pressure. The ATR absorbance using high pressure is about 10 times greater than with low pressure – all other sampling factors are identical. For rigid, crystalline, or hard, irregular surface samples we recommend a single reflection Diamond MIRacle ATR because it is easy to apply high pressure onto the small crystal (1.8 mm diameter) with the high-pressure clamp, producing over 10,000 psi.

The selection of the **ATR crystal characteristics** should be matched to the type of samples we run. Selection can be made to control depth of penetration of the IR beam, for hardness to prevent crystal damage, for desired spectral range and for acceptable pH range for acid or caustic samples. No individual crystal type will solve all problems, so PIKE Technologies offers a broad range of choices for ATR. Table 2 will give you some guidelines for selection of your ATR crystal.

**Table 2: ATR crystal characteristics for FTIR sampling.**

Crystal	n <sub>1</sub>	d <sub>p</sub> , for n <sub>2</sub> = 1.5 λ = 1000 cm <sup>-1</sup> , 45 deg, microns	Water Solubility g/100 g	pH Range	Hardness kg/mm
AMTIR	2.5	1.70	Insoluble	1–9	170
Diamond	2.4	2.01	Insoluble	1–14	5,700
Germanium	4.0	0.66	Insoluble	1–14	550
KRS-5	2.37	2.13	0.05	5–8	40
Silicon	3.4	0.85	Insoluble	1–12	1,150
ZnS	2.2	3.86	Insoluble	5–9	240
ZnSe	2.4	2.01	Insoluble	5–9	120

# PIKE Video Library Links —

## *Video Application Notes and Tips*

Click on the links below to access a wealth of informational videos featuring spectroscopic applications and sampling tips. Topics cover a range of sampling techniques including ATR, diffuse reflectance, transmission and more. Find other videos on our website such as PIKE software tutorials and an assortment of product showcases.



Proper Filling of an IR Liquid Cell



IR Transmission Measurements of Diffuse Reflectance




Accessing Formulation in Polymeric Materials



ATR: Single vs. Multiple Reflection



Troubleshooting Inverted Bands in the Spectrum

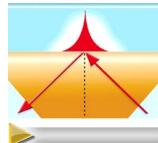

Depth of Penetration vs. Effective Penetration



Monitoring the Thermal Treatment of a Heterogeneous Catalyst



KBr: Proper Use and Handling Methods



Factors Influencing Spectral Quality for Mid-IR Diffuse Reflectance



The ABC's of Liquid Transmission Sampling



Calculating Critical Angle



Sample Preparation for Mid-IR Diffuse Reflectance